



149 Entin Road
Clifton, New Jersey 07014
www.nationalpublicseating.com

PSLT Series Assembly Instructions

Phenolic Science Lab table

Tools Required

WRENCH



Step 1: Place the tabletop upside-down on the floor with apron side up.

Note: Floor should be clear of any objects or dirt in order not to damage the tabletop.

Step 2: Remove plastic covers on bolts and place each leg in position with leg bolts going through the corner braces.

Step 3: Place a washer, lock washer and nut onto each bolt and tighten well using the provided wrench.

Step 4: Slide a leg boot onto each leg assuring that the wider side is facing the bottom of the leg.

Step 5: Screw a glide into the bottom of each leg.

Step 6: Turn the table over with the tabletop facing up. Place the table in desired location. Adjust the glides to assure that table is stable.

Step 7: Slide down protecting boots to cover the adjusting glides.

Dear Customer,

Thank you for your purchase. Although extreme care goes into manufacturing & packaging of this product, mistakes are possible. If you encounter any missing parts or difficulty in assembly, please contact our customer service Toll Free at 1-800-261-4112 or info@nationalpublicseating.com for courteous & immediate resolution to your problem.

CAUTION: Double check all screws to be tightened securely before use. Check all screws periodically.